

Name: Asma

Date: 09/30/2019

Goal:

1. Overnight cultures of pcb302 in A.Tume transformations from 9/26/19

Name: Asma

Date: 09/30/2019

Goal:

2. Overnight cultures of pcb302 in A.Tume transformations from 9/26/19

Protocol:

1. Add about 7 mL of YM to a 15 mL Falcon tube along with 7 μ L of kanamycin
2. Dip a p10 tip into your selected colony and drop into the tube
3. **Deviations:** Put 10ul of diH2O in eppendorf tube, picked a colony, and mixed colony with water very well. Added 2ul of this mix into falcon tube with 7ul kanamycin and 7ml YM.

*Placed cultures in shaking incubator @ 30 degrees @ 5pm on 09/30/2019 AK

*Colony+Water mix tubes labeled on long white tray in freezer

*Plates are parafilmed and refrigerated in new refrigerator

Conclusions:

Need to perform a colony pcr on samples.